making work easy

- DE
- EN
- FR
- IT
- ES
- PT
- TR
- PL
- RU
- ZH



# Mobiloskop S

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Made in Germany

21-6806 18112020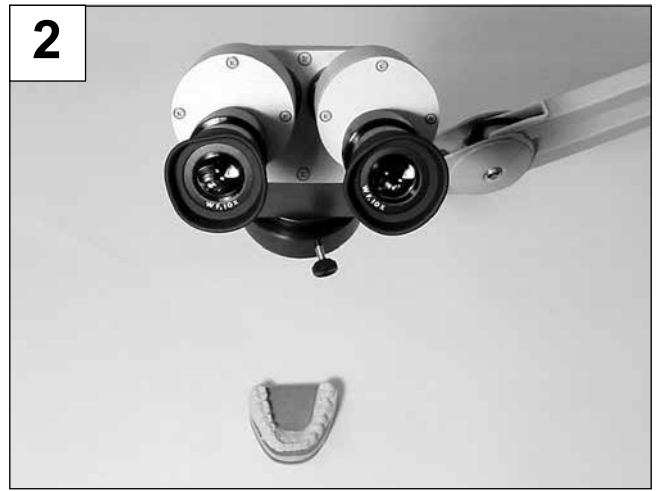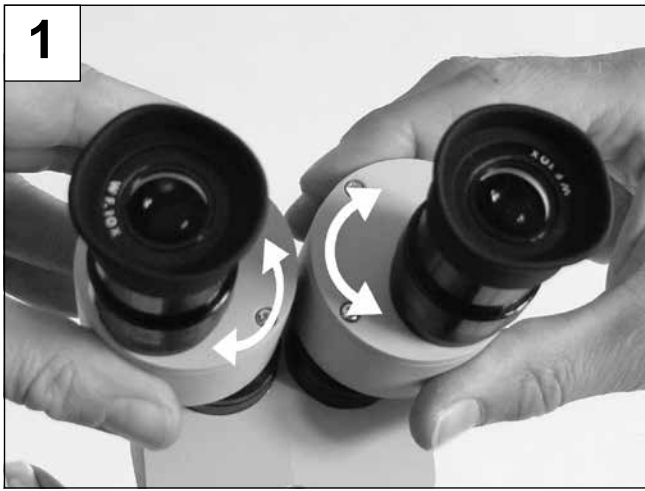

# Mobiloskop S

## Nr. 2200 0400

DEUTSCH

### 1. Einleitung

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ENGLISH

### 1. Introduction

We are pleased with your decision to purchase the Mobiloskop S, with swivel arm. The 5x or 10x magnification offers optimal detail control during precision tasks.

### 2. Application Area

The stereoscopic microscope with swivel arm is intended for work performed in laboratories, where it can be employed simultaneously at several workstations. The microscope's position is infinitely adjustable with the aid of the flexible swivel arm.

### 3. Hazards and warnings

- ▶ Remember that your field of view is greatly restricted when working with the Mobiloskop. Set up containers and nearby equipment (e.g., Bunsen burners) so that they cannot be knocked over.
- ▶ Working with magnifying equipment for long periods of time can impair your vision. Follow all guidelines issued by the responsible authorities (professional associations) and make sure you allow an adequate period of time for your eyes to readjust (judging distances)
- ▶ Information on REACH and SVHC is available on our website [www.renfert.com](http://www.renfert.com), in the Support area.

#### 3.1 Liability Exclusion

Renfert GmbH shall be absolved from all claims for damages or warranty if:

- ▶ The product is employed for any purposes other than those cited in the operating instructions;
- ▶ The product is altered in any way other than those alterations described in the operating instructions;
- ▶ The product is repaired by other than an authorized facility or if any but Renfert OEM parts are employed;
- ▶ The product continues to be employed, despite obvious safety faults.

### 4. Montage / Mise en service

#### 4.1 Installing the Device

For the installation of the support arm and of the microscope head, please refer to the installation instruction enclosed to the "Support arm with universal clamp".

#### 4.2 Adjusting the Focal Distance

Rotate the oculars to adjust the correct distance for your eyes (Figure 1).

#### 4.3 Adjusting the Dioptre Compensation

- ⇒ Place an object on the table top for orientation (Figure 2).
- ⇒ Turn the dioptre compensation sleeves completely clockwise (to their stops) (Figure 3).
- ⇒ Now turn the dioptre compensation sleeves one complete turn counter clockwise (Figure 4).
- ⇒ Place your left eye against the left eyepiece.
- ⇒ Use the left dioptre compensation sleeve to bring the object into sharp focus (Figure 5).
- ⇒ Place your right eye against the right eyepiece.
- ⇒ Use the right dioptre compensation sleeve to bring the object into sharp focus (Figure 6).

#### 4.4 Changing Magnification

Set the magnification to 5 x (0.5) or 10 x (1) (Figure 7).

### 5. Cleaning / Maintenance

Use a dry, lint-free cloth to clean the eyepiece and the protecting glass

**Never use solvents or cleansers containing abrasives!**

## 6. Spare Parts

You can find components subject to wear and the spare parts on the spare part list in the internet at [www.renfert.com/p918](http://www.renfert.com/p918).

The components excluded from the warranty (such as consumables or parts subject to wear and tear) are marked on the spare part list.

Serial number and date of manufacturing are shown on the type plate of the unit.

## 7. Guarantee

Provided the unit is properly used, Renfert warrants all components of the Mobiloskop S for a period of 3 years. The original sales receipt from the specialist dealer must be presented with any warranty claims. Components subject to natural wear are excluded from this warranty.

The warranty is voided in case of improper use; failure to observe the operating, cleaning, maintenance, and connection instructions; in case of independent repairs or repairs by unauthorized personnel; if spare parts from other manufacturers are employed, or; in case of unusual influences or influences not in compliance with the utilization instructions. Warranty service shall not extend the original warranty.

## 8. Technical Specifications

Light path:	45°, angled
Visual field diameter:	44 mm [1.734 inch] (5 x), 22 mm [0.867 inch] (10 x)
Focal depth:	22 mm [0.867 inch] (5 x), 10 mm [0.394 inch] (10 x)
Working distance, approx.:	150 mm [5.9 inch]
Weight, approx.:	1.8 kg [4.0 lbs]

## 9. Standard Delivery

- 1 Mobiloskop S
- 1 Support arm
- 1 Lens retainer
- 1 Lens with eyepieces
- 1 Protecting glass
- 1 Cover
- 1 Instruction manual

## 10. Delivery Versions

- No. 2200-0110 Eyepiece pair, 10 x
- No. 2200-0200 Additional 2 x lens, magnifies from 5 x / 10 x to 10 x / 20 x
- No. 3-0743 Replacement protecting glass to protect the lenses
- No. 2200-5000 Mobiloscope lighting
- No. 2200-0420 Stand with screw clamp

## 11. Error List

Problem	Possible cause	Corrective action
<b>No visible image.</b>	<ul style="list-style-type: none"> <li>• Magnification change knob incorrectly adjusted.</li> </ul>	<ul style="list-style-type: none"> <li>• Turn the knob to its end stop.</li> </ul>
<b>Image is not sharply focussed.</b>	<ul style="list-style-type: none"> <li>• Incorrect dioptre compensation setting.</li> <li>• Incorrect distance to the object.</li> <li>• Dirty lens / protecting glass.</li> </ul>	<ul style="list-style-type: none"> <li>• Repeat the dioptre compensation procedure.</li> <li>• Change the distance to the object.</li> <li>• Use a clean, lint free cloth to clean the lens / protecting glass.</li> </ul>
<b>The support arm does not stand still.</b>	<ul style="list-style-type: none"> <li>• Fixation is loose.</li> </ul>	<ul style="list-style-type: none"> <li>• Tighten fixation.</li> </ul>

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Renfert GmbH • Untere Gießwiesen 2 • 78247 Hilzingen/Germany