making work easy

- FAQ - Deutsch
- FAQ - English
- FAQ - Français
- FAQ - Italiano
- FAQ - Italiano
- FAQ - Português
- FAQ - Türkçe
- FAQ - Русский
- FAQ - Polski
- FAQ - Čeština
- FAQ - Українська
- FAQ - 中文
- FAQ - 日本語
- FAQ - 한국어
- FAQ - اللغة العربية
- FAQ - Svenska

# Vibrax

FAQ - frequently asked questions

Made in Germany

# FAQ - Deutsch

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# FAQ - English

*For chapter indications, refer to the complete instruction manual on the internet.*

Error	Possible cause	Corrective action
<p><b>The device generates no vibrations.</b></p>	<ul style="list-style-type: none"> <li>• Power cord not properly plugged in at the device.</li> <li>• Blown fuse.</li> <li>• No power at the wall outlet.</li> </ul>	<ul style="list-style-type: none"> <li>• Check that the power cord is securely seated in the socket.</li> <li>• Replace the fuse (refer to Sec. 3.4).</li> <li>• Test the wall outlet for power.</li> </ul>
<p><b>No difference in the vibration intensity between the switch settings Fine / Coarse.</b></p>	<ul style="list-style-type: none"> <li>• Faulty diode (short circuit).</li> </ul>	<ul style="list-style-type: none"> <li>• Return the device for service.</li> </ul>
<p><b>No function in coarse oscillation mode.</b></p>	<ul style="list-style-type: none"> <li>• Faulty diode (interruption).</li> <li>• Faulty switch.</li> <li>• Faulty switch connection wiring.</li> </ul>	<ul style="list-style-type: none"> <li>• Return the device for service.</li> </ul>
<p><b>No function in fine oscillation mode.</b></p>	<ul style="list-style-type: none"> <li>• Faulty switch.</li> <li>• Faulty switch connection wiring.</li> </ul>	<ul style="list-style-type: none"> <li>• Return the device for service.</li> </ul>
<p><b>No effect at the different power level settings.</b></p>	<ul style="list-style-type: none"> <li>• Faulty switch.</li> <li>• Faulty switch connection wiring.</li> </ul>	<ul style="list-style-type: none"> <li>• Return the device for service.</li> </ul>
<p><b>Device extremely loud during operation.</b></p>	<ul style="list-style-type: none"> <li>• Loose internal components.</li> <li>• Loose rubber mat.</li> </ul>	<ul style="list-style-type: none"> <li>• Return the device for service.</li> <li>• Attach the rubber mat correctly (see Chap. 1.2).</li> </ul>

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